

## **FORMALIN VAPORISER**



The CAS Formalin Vaporiser is designed as a compact and simple unit to enable safety cabinets to be fumigated using formaldehyde.

The unit can be fitted as an integral part of the cabinet or alternatively supplied as a free standing unit that is placed inside the cabinet and plugged into an internal socket. If fitted to a CAS Safety Cabinet the unit can be wired independently or directly into the cabinet controls to allow the vaporiser to boil off the fumigant over a timed period.

To run the unit, remove the large cap on the top and pour in the required quantity of formalin with an equal quantity of water; replace the cap, then set the unit in operation to boil off the fumigant.

## **Technical Data**

Size: 185mm wide x 105mm deep x 185mm high

Capacity: 60ml

Electrical Supply: 230v/1ph/50Hz

Electrical Load: 2 amps